
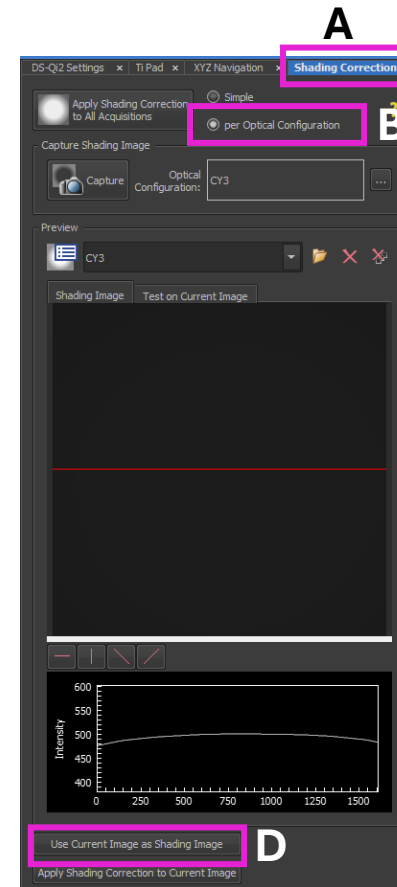


Set up shading correction for fluorescence imaging in Elements

1. Open shading correction image
 - Go to Shading Correction > General Use folder in D:\ drive
 - Choose the objective lens you are using (4X, 10X, 20X)
 - Select shading correction image based on background pixel number and fluorescence channel
2. Go to Shading correction tab (A)
3. Make sure “per optical configuration” is selected (B)
4. Choose the fluorescent channel you need for the shading correction by clicking the ‘...’ (C)
5. Click “Use Current Image as Shading Image” (D)
6. **Make sure you turn on shading correction by clicking the button left of the ‘play’ button (E)**

Note: Only need to do this once at the beginning of imaging session!

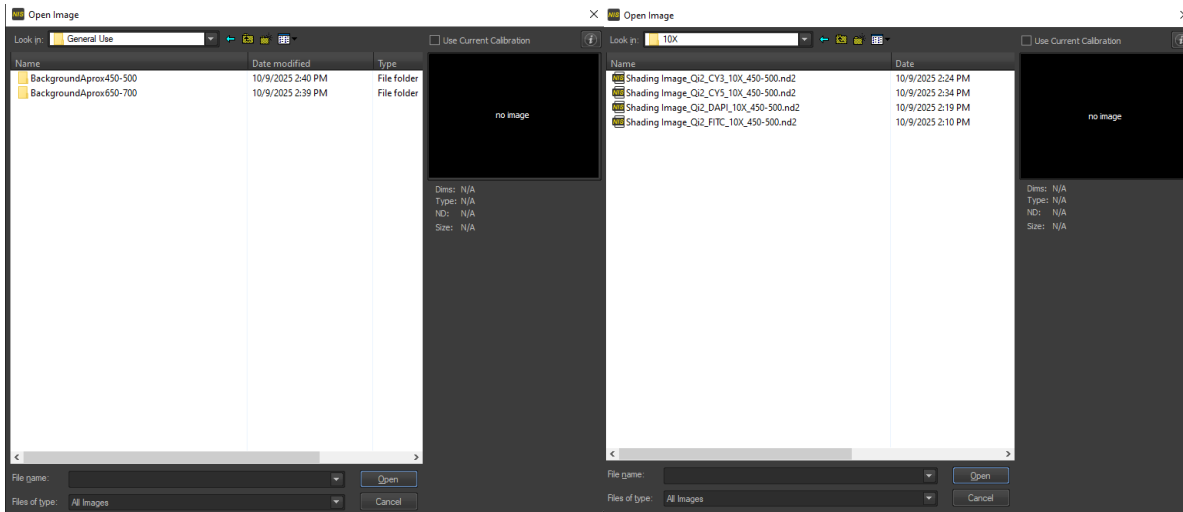

Off:



On:



A

B

C

D

E